

# "The Sealed, Safe & Efficient Way" *Recommended for Mice in Transit*

Wouldn't it be nice if shipping containers were:

- ☐ divided into sealed compartments so that species and sexes couldn't intermingle
- ☐ easy to remove animals avoiding bites and scratches
- ☐ designed to avoid stress and injury to the animal
- ☐ had lots of view ports and air holes to avoid unnecessary opening
- ☐ came sterile or *on site* autoclaveable and could be directly filled in a barrier
- ☐ could be easily unpacked (no handling) with mice sliding right into their permanent cages



We Thought so too; and also added: **100% recyclable!**


## The **TRANS-TAINER** system: **Tube, Caps, Carrier & Hydration Gel**

### THE TUBE

114 grams live animal weight

The shape rodents feel comfortable in, they build tunnels! No corners or any place to begin gnawing. Reusable Opaque Polypropylene & Clear Polycarbonate are easily sanitized in cage washers and are safe for temperatures and disinfectants found in most facilities.

Simple, cost effective and economical! Barrier Friendly - Entry into the barrier is a simple spray or soak.

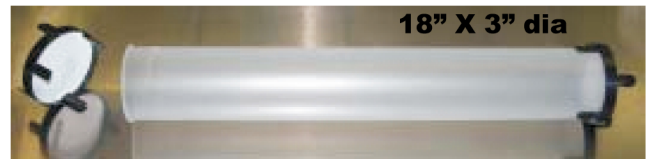

Available in 9" tube for "short term" inter facility travel.



### THE END CAP

Securely snaps on, seal must be broken to remove. When cap is removed tube is slightly tilted and animals slide gently into cage. A new cap is always used and the .5 Micron polypropylene filter is ideal for hazardous species and reduces risk to handlers

### THE CARRIER BOX

The Box will take three separate Trans-Tainer tubes effectively separating strains, species, micro-flora, health status & sexes. Tube isolation prevents cross contamination. Assembly and loading instructions are clearly printed on box. "LIVE ANIMAL" alerts are very visible to all courriers and handlers. Flaps fold down to create tabs which create an "air gap" so that end of box cannot be blocked off.

22.5" X 17.5" X 7.



**TRANS-TAINER®**

**SPECIALTY MOUSE SHIPPER**

[www.perotech.com/transainer/](http://www.perotech.com/transainer/)

IN CANADA:

**Perotech Sciences Inc**


55 Albert St., Markham ON L3P 2T4

[www.perotech.com](http://www.perotech.com)

416-332-8394


[info@perotech.com](mailto:info@perotech.com)

# The Convenient Reliable Moisture Source *Recommended for Mice in Transit*

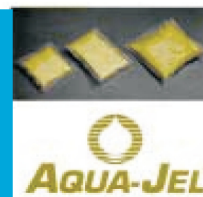

## AQUA-JEL

The gel is a reliable moisture source to prevent dehydration. It will not spill or get the animals wet - an amazing 98% available to the animal. Aqua-Jel is commercially sterile until opened and can be further gamma irradiated. No refrigeration is required, it's easy to use and store and comes in 3 convenient sizes.

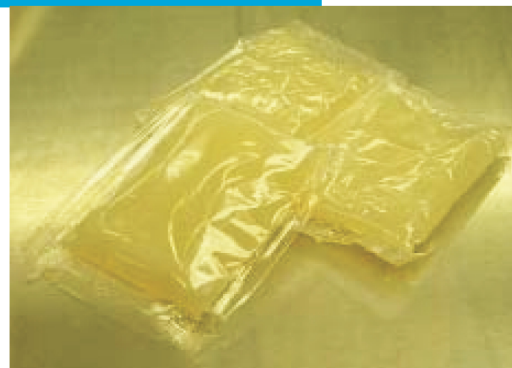

### ORDERING INFORMATION

MODEL	SIZE	WEIGHT	PACKS/CASE
86-10	86 gram	1.5 lb	10
86-50	86 gram	7 lb	50
86-150	86 gram	21lb	150
150-100	150 gram	23.5 lb	100
300-50	300 gram	47 lb	50

## TRANS-TAINER M10-3 ORDERING INFORMATION

MODEL	DESCRIPTION & SIZE	PACKS/CASE
CT-3	Corrugated Carrier Knocked Down complete with Instructions, Will Hold 3 Tubes	10
PPC	Opaque Polypropylene Tubes With Filtered End Caps	10
PCC	Clear Polycarbonate Tubes With Filtered End Caps	10
TC-16	Polysulphone Transfer Tube with Filtered End Caps	10
EC-G	End Caps with Filter, Packaged 6 Per Autoclave Bag	6
M10-PS Kit	Complete with Carrier, 3 Tubes, 3 86g Aqua-Jel pouches	1

\* **Phone For:** Single Case, US Dollar, FOB FACTORY, VOLUME, BLANKET ORDER AND CONTRACT PRICING



< Calculate how long I'll be in transport and then cap me in with Jel, food & bedding. At my destination after tube is received to protocol, I'm taken to my cage. One end cap is easily snapped open and the tube is tilted and I slide safely into my new cage. >

**416 332-8394**

